

Product Note M165 - 01/17

B-KIMW ATR-IR Polymers, Plastics and Additives Library, Kunststoff-Institut Lüdenscheid

With a production volume of hundreds of megatons each year, polymers and plastics have become one of the most important basic materials for manufacturing. To always obtain the required material properties many different basic polymer types are available and new types are constantly developed. Moreover, by blending basic polymers and adding fillers and additives an almost countless number of individual plastic materials is created.

The key to effective polymer identification

Out of the various analytical methods used for the investigation of polymers and polymer products, FTIR (Fourier transform infrared) spectroscopy is among the fastest and most powerful tools available. Besides basic polymers, also fillers, additives and other components contribute to the measured IR sample spectrum. Hence, the IR spectrum provides a chemical fingerprint of a certain plastic material.

With the availability of reference spectra databases, identification of plastic materials can be performed within just a minute. The most efficient and convenient method to obtain an IR spectrum is the attenuated total reflection (ATR) technique since generally no or only little sample preparation is needed and results are available in seconds.



Certified Samples. Quality Data. Assembled by Experts

The "Polymers, Plastics and Additives Library" of the Kunststoff-Institut Lüdenscheid is an extensive collection of polymer ATR spectra.



A close collaboration between the Kunststoff-Institut Lüdenscheid (KIMW) and Bruker led to high quality spectra of currently used plastic materials including technical biopolymers. All materials were carefully verified and selected by the KIMW and are kept as retained reference samples.

First Quality, then quantity

No other spectral library offers such quality standards and high information content for each entry. Each sample is well-characterized by its plastic type and trade name. Furthermore, the library contains a wealth of additional knowledge suchs as manufacturer, color, fillers, application area and physical properties.

Polymer PBT CAS Registry Number 26062-94-2 Substance polybutylene terephthalate Trade Name Celanex 2002 Supplier Celanese Filler unfilled Filler Content 0 Color natural Processing Method injection molding Metting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	Compound informati	on			
CAS Registry Number 26062-94-2 Substance polybutylene terephthalate Trade Name Celanex 2002 Supplier Celanese Filler unfilled Filler Content 0 Color natural Processing Method injection molding Metting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunstsoff Institut Lüdenscheid 0140	1. A.		PBT		
Substance polybutylene terephthalate Trade Name Celanex 2002 Supplier Celanese Filler unfilled Filler Content 0 Color natural Processing Method injection molding Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140		er			
Trade Name Celanex 2002 Supplier Celanese Filler unfilled Filler Content 0 Color natural Processing Method injection molding Metting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	and the second se			alate	
Supplier Celanese Filler unfilled Filler Content 0 Color natural Processing Method injection molding Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	STREET STREET				
Filler unfilled Filler Content 0 Color natural Processing Method injection molding Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Reference Kunststoff Institut Lüdenscheid 0140					
Color natural Processing Method injection molding Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140					
Color natural Processing Method injection molding Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	and the second		110		
Processing Method injection molding Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140			•		
Melting Point 225 Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140					
Glass Transition 60 Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	and the second				
Young's modulus (RT) 2600 MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140					
MVR/MFR 20 cm3/ 10 min [2,16 kg/260 °C] Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140		T)			
Density 1.31 Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	and the second	.,		kg/260 °C1	
Hardness (Shore) 78 [Rockwell M] Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	Density		and the second sec		
Measurement Technique DiamondATR Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140	a hand a hand a second s		78 [Rockwell M]		
Sample Preparation cut surface, core material Sample Reference Kunststoff Institut Lüdenscheid 0140		ique			
Sample Reference Kunststoff Institut Lüdenscheid 0140			cut surface, core mat	erial	
			Kunststoff Institut Lüd	enscheid 0140	
				nı l	
	00- 		2000 1500 ality Compound Name PBT	1000	
2 979 PBT 26062-94-2	4000 3000 Hit No. ▲ Color	Hit Qua	ality Compound Name PBT	CAS number 26062-94-2	

Figure 1: Result of spectrum identification with the ATR-IR Polymers, Plastics and Additives Library. The upper window shows extensive compound information as well as how the reference spectra were collected.

Some of the supplementary information provided by the KIMW-library is shown in the upper part of Figure 1. To always provide a representative spectrum of each material several spectra were measured at different sample positions and averaged for the final library entry. In cases were the surface material differed from the core, both spectra were added to the database.

To guarantee highest quality spectra, measurements were done on a Bruker TENSOR II spectrometer equipped with a platinum diamond ATR and a spectral range down to 340 cm⁻¹. Furthermore, dark samples were additionally measured with a germanium ATR unit to suppress artifacts resulting from fillers with a high refractive index like carbon black. The database is completed by spectra of relevant polymer additives like fillers, stabilizers, plasticizers, antistatic agents etc.

Always Up-to-Date

As new plastic materials are entering the market also reference databases have to be updated regularly. Therefore, the "Polymers, Plastics and Additives Library" will be continuously expanded by new materials. In case of specific questions regarding samples and spectra which are included in the library, the Kunststoff-Institut Lüdenscheid provides support.

Kunststoff-Institut Lüdenscheid (KIMW)

- Leading provider of in-depth services in a wide range of polymer technologies
- Supports customers in selecting, developing, optimizing and implementing products, tools and processes in all areas of plastics technology.
- DIN EN ISO 9001 certified
- DIN EN ISO/IEC 17025:2000 accredited laboratory
- Internet: www.kunststoff-institut.de

Overview:

B-KIMW

Bruker ATR-IR Polymers, Plastics and Additives Library, Kunststoff-Institut Lüdenscheid

Specifications:

- Around 1050 high quality ATR spectra of up-to-date plastic materials including technical biopolymers, and additives
- Sample materials selected and analyzed by the polymer institute KIMW
- Ca. 160 different polymer types and about 150 additives
- Spectral quality and reference information verified by the polymer institute KIMW
- Comprehensive material information from material data sheets
- Full MIR range: 4000 340 cm⁻¹
- Additional Germanium-ATR spectra of materials with high refractive index
- Regular library update available with additional new plastic materials and additives (once per year).

www.bruker.com/optics

Bruker Optics

Bruker Optics is continually improving its products and reserves the right to change specifications without notice. © 2018 Bruker Optics BOPT - 4001031 - 02